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## Instruction for Use TPP TubeSpin® Bioreactor

The TubeSpin® bioreactor is suitable for high-throughput screening, cultivation and optimization processes of prokaryotic (aerobic bacteria) and eukaryotic (yeasts, algae, HEK, CHO, Sf-9) suspension cells.

The TubeSpin® bioreactor is available with a filter screw cap with 5 or 10 openings over a PTFE filter membrane.

The TubeSpin® Bioreactor is intended for single use only.

### Note:

Follow national regulations when handling biological material, use appropriate protective clothing


Take note of the regulations of aseptic working during the work process.

### Instructions:

- Check the packaging material and the product for intactness. Use only faultless material.
- Check the expiry date (EXP) on label and packaging. Only use products with valid EXP.
- Open the packaging in a sterile environment and remove an entire system for use.
- Open the bioreactor and fill it with the medium and inoculum according to your laboratory routine. Please refer to the optimal filling quantity (see technical data).
- Close the filter screw cap with ventilation holes. A continuous gas exchange takes place through the integrated hydrophobic membrane 0.22  $\mu\text{m}$ .

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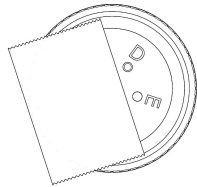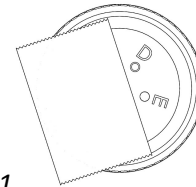
- *If the aeration/gas exchange needs to be reduced, individual openings may be closed using lab appropriate labels or tape (picture 1)*

• Picture 1

- **Cultivate:** Shake cells in appropriate incubation shakers. Refer to the cultivation conditions of the cells (temperature, CO<sub>2</sub> and humidity control, shaking frequency)
- **Shakers:** Follow the safety instructions and operating instructions of the shaker manufacturer.
- **Centrifuging:** The use of suitable rotors or centrifuge adapters is recommended. Follow the safety instructions of the centrifuge manufacturer.
- **The mechanical strength of the tubes during centrifugation is influenced by**
  - Shape and material
  - Accuracy of fit in the centrifuge adapter
  - Temperature, centrifugation time, g-number
  - Chemical and physical properties of the centrifuged material
  - Rotor type: fixed-angle or swing-out rotor
- **Test the multitude of influencing factors under routine conditions beforehand.**
- **Storage before use at room temperature, protected from UV light**

#### Remarks and Recommendations:

- *The graduation serves as a guide for filling quantity and is no absolute value.*
- *Alcohol 90 % with mechanical stress (rubbing) might dissolve or blur the white color of the print.*

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#### Additional:

Instructions for use, chemical resistance lists, and quality certificates of the respective products are available for download on the homepage [www.tpp.ch](http://www.tpp.ch).

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